Scan QR code to download  
the full instruction manual.

# Adventurer™ Balances User Guide

## REVISION HISTORY

| Date               | Version | Description   |
|--------------------|---------|---|
| March 9, 2023      | B       | We incorporate the product specification table into the instruction manual in an effort to reduce printed materials.<br>Please download the user instruction manual at <a href="https://ohaus.com/products">ohaus.com/products</a> or use your smartphone to scan the QR code on the front page of this manual. |
| September 27, 2023 | C       | Updated company address   |

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## 1. INTRODUCTION

Thank you for purchasing our product. This user guide will help you install your balance and use it safely. Please read it carefully before installation and operation. For detailed instructions, please visit [www.ohaus.com/products](http://www.ohaus.com/products).

## 2. SAFETY INFORMATION

This manual contains installation, operation and maintenance instructions for AX Series balance. Please read this manual completely before installation and operation.

### Definition of Signal Warnings and Symbols

Safety notes are marked with signal words and warning symbols. These show safety issues and warnings. Ignoring the safety notes may lead to personal injury, damage to the instrument, malfunctions and false results.

**WARNING** For a hazardous situation with medium risk, possibly resulting in injuries or death if not avoided.

**CAUTION** For a hazardous situation with low risk, resulting in damage to the device or the property or in loss of data, or injuries if not avoided.

**Attention** For important information about the product

**Note** For useful information about the product

Warning Symbols

General Hazard

Electrical Shock Hazard

Explosion hazard

### Safety Precautions

**CAUTION:** Read all safety warnings before installing, making connections, or servicing this equipment. Failure to comply with these warnings could result in personal injury and/or property damage. Retain all instructions for future reference.

- Before connecting power, verify that the AC adapter's input voltage range and plug type are compatible with the local AC mains power supply.
- Do not position the equipment such that it is difficult to reach the power connection.
- Make sure that the power cord does not pose a potential obstacle or tripping hazard.
- Operate the equipment only under ambient conditions specified in these instructions.
- This equipment is for indoor use only.
- Do not operate the equipment in wet, hazardous or unstable environments.
- Do not allow liquids to enter the equipment.
- Do not load the equipment above its rated capacity.
- Do not drop loads on the platform.
- Do not place the equipment upside down on the platform.
- Use only approved accessories and peripherals.
- Disconnect the equipment from the power supply when cleaning.
- Service should only be performed by authorized personnel.

**WARNING:** Electrical shock hazards exist within the housing. The housing should only be opened by authorized and qualified personnel. Remove all power connections to the unit before opening.

**WARNING:** Never work in an environment subject to explosion hazards! The housing of the instrument is not gas tight. (Explosion hazard due to spark formation, corrosion caused by the ingress of gases).

### Intended Use

This instrument is intended for use in laboratories, pharmacies, schools, businesses and light industry. It must only be used for measuring the parameters described in these operating instructions. Any other type of use and operation beyond the limits of technical specifications, without written consent from OHAUS, is considered as not intended.

This instrument complies with current industry standards and the recognized safety regulations; however, it can constitute a hazard in use.

If the instrument is not used according to these operating instructions, the intended protection provided by the instrument may be impaired.

## 3. INSTALLATION

### 3.1 Select the Location

The location must be sturdy, flat and level.

Avoid locations with excessive air current, vibrations, heat sources or rapid temperature changes.

Allow sufficient space around the instrument.

### 3.2 Leveling the Equipment

To level the instrument, adjust the feet/leveling wheel so the bubble is centered in the circle of the level indicator. Be sure the equipment is level each time its location is changed.

### 3.3 Connecting Power

Connect the AC adapter power cord to the instrument's power input connector, and then connect the AC plug to a suitable electrical outlet.

**Attention:** Only use an AC adapter specified by OHAUS.

**Attention:** For optimal weighing performance, allow the balance to warm up for 60 minutes prior to use.

### 3.4 Connecting the Interface

Use the built-in RS-232 Port to connect either to a computer or a printer with a standard (straight-through) serial cable. Or connect using the scale's USB port.

Interface connections on the rear of the balance:

USB

RS232

USB1: Used to connect to PC only

USB2: Used to connect to a USB flash driver only

RS232: Used to connect to PC or Printer

USB connection on the front of the balance:

### 3.5 Initial Calibration

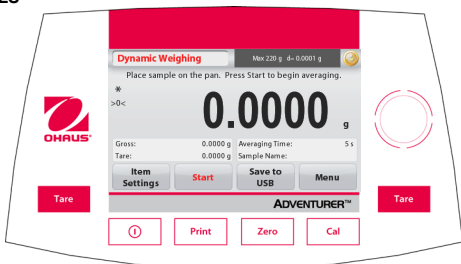
For best results, the instrument's calibration should be checked prior to first use. If adjustment is needed, refer to the Calibration section of the instruction manual.

## 4. OPERATION

### 4.1 Overview of Display, Home Screen

This equipment utilizes a touch-sensitive display. Touch areas and Buttons to control the equipment's functions.

#### CONTROLS

| Button   | Action  |
|--|---|
|  | User logout button  |
|  | Short Press (if powered Off): Turns on the scale<br>Long Press (if powered On): Turns off the scale<br>Note: The balance will automatically power on when power is connected. |
|  | Short Press: Prints the present data to a printer or a computer.  |
|  | Short Press: Perform Zero operation   |
|  | Short Press: Perform Calibration operation  |
|  | Short Press: Perform Tare operation   |

## Main Application Screen

Application

Instructional Messages  
Stability (\*), Net (NET),  
Gross (G) and/or center  
of zero (>0<) indicators

Reference Fields

User logout button

Result Field: Information varies by  
application  
Touch g to change unit

Application Buttons:  
Functions vary by application

## MENU & SCREEN NAVIGATION

Touch **Menu** to open the menu list.

### Calibration:

Touch to view calibration options.

### Balance Setup:

Touch to view and change balance settings.

### Weighing Units:

Touch to view and change weighing units.

### Data Maintenance:

Touch to view data maintenance settings.

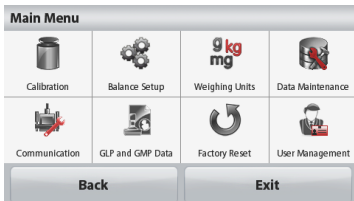
### Communication:

Touch to view COM Device Settings and Print Settings.

### GLP and GMP Data:

Insert user data for traceability.

### Factory Reset:

Touch to do a Factory reset of menu settings.

### User Management:

Touch to enter the submenus:  
**User Profiles** (Edit the user),  
**Change Password** and **Auto Standby**

## 4.2 Using the Balance

Note: Before using any application, be sure the balance has been leveled and calibrated.

### Weighing Application

1. If required, place an empty container on the pan and press Tare.
2. Add sample to the pan or container. The display shows the weight of the sample.

## 5. MAINTENANCE

### 5.1 Cleaning

**WARNING:** Electric Shock Hazard. Disconnect the equipment from the power supply before cleaning.

Make sure that no liquid enters the interior of the instrument.

**Attention:** Do not use solvents, harsh chemicals, ammonia or abrasive cleaning agents.

Glass surfaces may be cleaned with a commercial glass cleaner.

The housing may be cleaned with a damp cloth and mild detergent if necessary.

## 5.2 Service Information

For troubleshooting common problems, download the user manual from [www.Ohaus.com](http://www.Ohaus.com). If the issue persists, please visit our website for available support and service centers at <https://www.ohaus.com/en-us/support/service-centers>.

## 6. TECHNICAL DATA

The technical data is valid under the following conditions:

- Indoor use only
- Altitude: Up to 2000 m
- Operating temperature: 5°C to 40°C
- Operating temperature for optimum weighing performance: 10°C to 30°C
- Humidity: maximum relative humidity 80% for temperatures up to 31°C, decreasing linearly to 50% relative humidity at 40°C
- Electrical supply: 12VDC, 0.84A. (For use with certified or approved power supply, which must have a SELV and limited energy output.)
- Mains supply voltage fluctuations: up to  $\pm 10\%$  of the nominal voltage
- Installation category II
- Pollution degree: 2

Detailed product specifications can be found in the user manual and can be downloaded from [www.Ohaus.com/products](http://www.Ohaus.com/products).

### LIMITED WARRANTY

Ohaus products are warranted against defects in materials and workmanship from the date of delivery through the duration of the warranty period. During the warranty period Ohaus will repair, or replace any component(s) that proves to be defective at no charge, provided that the product is returned, freight prepaid, to Ohaus.

This warranty does not apply if the product has been damaged by accident or misuse, exposed to radioactive or corrosive materials, has foreign material penetrating to the inside of the product, or as a result of service or modification by other than Ohaus. In lieu of a properly returned warranty registration card, the warranty period shall begin on the date of shipment to the authorized dealer. No other express or implied warranty is given by Ohaus Corporation. Ohaus Corporation shall not be liable for any consequential damages.

As warranty legislation differs from state to state and country to country, please contact Ohaus or your local Ohaus dealer for further details.

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P/N 30283442 C © 2023 Ohaus Corporation, all rights reserved.