Scan QR code to download
the full instruction manual.

PJX Series Balances User Guide

REVISION HISTORY

Date	Version	Description
November 25,2022	F	We incorporate the product specification table into the instruction manual in an effort to reduce printed materials. Please download the user instruction manual at ohaus.com/products or use your smartphone to scan the QR code on the front page of this manual.
September 12, 2023	G	Updated the company address

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1. INTRODUCTION

Thank you for purchasing our product. This user guide will help you install your balance and use it safely. Please read it carefully before installation and operation. For detailed instructions, please visit www.ohaus.com/products.

2. SAFETY INFORMATION

This manual contains installation, operation and maintenance instructions for PJX Series balance. Please read this manual completely before installation and operation.

Definition of Signal Warnings and Symbols

Safety notes are marked with signal words and warning symbols. These show safety issues and warnings. Ignoring the safety notes may lead to personal injury, damage to the instrument, malfunctions and false results.

WARNING For a hazardous situation with medium risk, possibly resulting in injuries or death if not avoided.

CAUTION For a hazardous situation with low risk, resulting in damage to the device or the property or in loss of data, or injuries if not avoided.

Attention For important information about the product

Note For useful information about the product

Warning Symbols

General Hazard

Electrical Shock Hazard

Alternating current

Direct current

Safety Precautions

CAUTION: Read all safety warnings before installing, making connections, or servicing this equipment. Failure to comply with these warnings could result in personal injury and/or property damage. Retain all instructions for future reference.

- Verify that the AC adapter's input voltage range and plug type are compatible with the local AC mains power supply.
- Do not position the equipment such that it is difficult to reach the power connection.
- Make sure that the power cord does not pose a potential obstacle or tripping hazard.
- Operate the equipment only under ambient conditions specified in these instructions.
- This equipment is for indoor use only. Do not operate the equipment in hazardous or unstable environments.
- Do not drop loads on the platform.
- Use the balance only in dry locations.
- Use only approved accessories and peripherals.
- Disconnect the equipment from the power supply when cleaning.
- Service should only be performed by authorized personnel.

Intended Use

Use the instrument exclusively for weighing as described in the operating instructions. Any other type of use and operation beyond the limits of technical specifications without written consent from OHAUS, is considered as not intended.

This instrument complies with current industry standards and the recognized safety regulations; however, it can constitute a hazard in use.

If the instrument is not used according to these operating instructions, the intended protection of the instrument may be compromised and OHAUS assumes no liability.

3. INSTALLATION


3.1 Select the Location


Avoid heat sources

Rapid temperature changes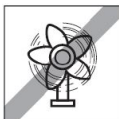

Air current or excessive vibrations

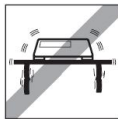Allow sufficient space


3.2 Leveling the Equipment

Be sure the balance is level before it is used or after its location is changed.

The PJX balance has a level bubble in a small round window beside the display.

To level the balance, adjust the 4 leveling feet until the bubble is centered in the circle.

Please refer to the right figure for leveling.

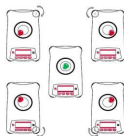3.3 Connecting Power and Acclimatising the Balance

Connect the AC adapter power cord to the instrument's power input connector, and then connect the AC plug to a suitable electrical outlet.

Acclimatising

It is suggested that the balance should not be used until it has been connected to power and acclimatised to the environment for a certain period of time. In the case of a balance with the precision above 0.1 mg, the acclimatisation time should be 1.5 hours; in the case of balance with the precision of 0.01 mg, the acclimatisation time should be more than 4 hours.

3.4 Connecting the Interface

The PJX balance has two connectivities, RS232 port and Mini DIN.

Use the RS-232 port to connect either to a computer or a printer with a standard (straight-through) serial cable.

Mini DIN is used to connect with another accessory, as an alternative option for RS232. Interface connections on the rear of the balance.

Mini DIN

RS232

RS232: Used to connect to PC or Printer

Note: See the Printing section for Connecting, Configuring and Testing the Printer / Computer Interface.


3.5 Initial Calibration

For best results, the instrument's calibration should be checked prior to first use. If adjustment is needed, refer to the Calibration section of the instruction manual.

4. OPERATION

4.1 Overview of Display, Home Screen Controls

Controls

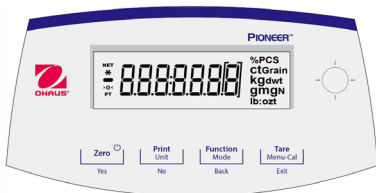CONTROL FUNCTIONS

Button	Zero	Print Unit	Function Mode	Tare Menu-Cal
	Yes	No	Back	Exit
Primary Function (Short Press) 	On / Zero - If the balance is Off, turns on the balance. - If the balance is On, sets zero.	Print Sends the current displayed value to the serial interface.	Function Operation is dependent on the application mode.	Tare Performs tare operation.
Secondary Function (Press and Hold) (Extended Press)	Off - Zeros the current value. - None	Unit - Changes weighing units. - None	Mode - Changes application mode. - None	Menu-Cal - Enters the main menu. Calibration is the first sub-menu. - Views the preset Tare value.
Menu Function (Short Press) 	Yes Accepts the current (blinking) setting on the display.	No - Rejects the current (blinking) setting on the display. - Increments a value being entered.	Back - Reverts back to previous menu item. - Decrements a value being entered.	Exit - Immediately exits the sub-menu. - Aborts a calibration in progress.

Display

Net (NET)
Stability (*)
Negative (-)
Centre of Zero (>0<)
Preset tare (PT)

Result Field: depending on application

Unit

4.2 Principal Function

Zeroing: Press **Zero** to zero the balance.

Taring: With no load on the pan, press **Zero** to set the display to zero. Place an empty container on the pan and press **Tare**. Add objects to the container and its net weight is displayed. After the container and the objects are removed, the tare value will be displayed as a negative number. Press **Tare** to clear.

Weighing: First press **Zero** to set the display to zero. Place an object on the pan. The display indicates the gross weight.

5. MAINTENANCE

5.1 Cleaning

WARNING: Disconnect the equipment from the power supply before cleaning.

Make sure that no liquid enters the interior of the balance.

Clean the Balance at regular intervals.

Housing surfaces may be cleaned with a lint-free cloth slightly dampened with water or a mild cleaning agent.

Glass surfaces may be cleaned with a commercial glass cleaner.

Attention: Do not use solvents, harsh chemicals, ammonia or abrasive cleaning agents.

5.2 Service Information

For troubleshooting common problems, download the user manual from www.Ohaus.com. If the issue persists, please visit our website for available support and service centers at <https://www.ohaus.com/en-us/support/service-centers>.

6. TECHNICAL DATA

The technical data is valid under the following conditions:

- Indoor use only
- Altitude: Up to 2000 m
- Operating temperature range: 5°C to 40°C
- Operating temperature for optimum weighing performance: 10°C to 30°C
- Humidity: maximum relative humidity 80% for temperatures up to 30°C, decreasing linearly to 50% relative humidity at 40°C
- Mains supply voltage fluctuations: up to $\pm 10\%$ of the nominal voltage
- Installation category II
- Pollution degree: 2

Detailed product specifications can be found in the user manual and can be downloaded from www.Ohaus.com/products.

LIMITED WARRANTY

Ohaus products are warranted against defects in materials and workmanship from the date of delivery through the duration of the warranty period. During the warranty period Ohaus will repair, or replace any component(s) that proves to be defective at no charge, provided that the product is returned, freight prepaid, to Ohaus.

This warranty does not apply if the product has been damaged by accident or misuse, exposed to radioactive or corrosive materials, has foreign material penetrating to the inside of the product, or as a result of service or modification by other than Ohaus. In lieu of a properly returned warranty registration card, the warranty period shall begin on the date of shipment to the authorized dealer. No other express or implied warranty is given by Ohaus Corporation. Ohaus Corporation shall not be liable for any consequential damages.

As warranty legislation differs from state to state and country to country, please contact Ohaus or your local Ohaus dealer for further details.

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P/N 30431477 G © 2023 Ohaus Corporation, all rights reserved.